

Instructions for GazFit Connectors

Make sure you have the correct sized GazFit for your torch. GazFit connectors are available to suit 8mm, 6mm, 5mm and 4mm side arms.

Slide the flexible argon tubing over the barbed end of the GazFit connector.

Slide the end with the o-rings over the torch side arm.

NOTE: Since the GazFit connector expands slightly to fit the torch side arm and there is a small variation in the diameter of the side arms, the GazFit connector may appear loose when it is moved from one torch to another. If this is the case, let the GazFit connector rest for 24 hours and it will relax to its original diameter. This process can be speeded up by putting the GazFit connector in hot water. If the connection is still loose, replace the GazFit connector with a new one.