



GLASS EXPANSION

Quality By Design

Instructions for Nexus Universal Nebulizer Connection Kit

Part Number FT-16-8-X


General Safety

- It is recommended you read and understand this procedure before starting.
- It is advised you wear safety gloves and glasses when handling glass.


Disclaimer: Glass Expansion supplies products to customers with the assumption that only qualified individuals who are properly trained in the appropriate procedures will use our products. It is our customer's entire responsibility to safely store, handle and use the products ordered from Glass Expansion. We assume the requisitioner to be experienced, competent, safety conscious professionals. Glass Expansion assumes no liability for any damage or injuries -physical, emotional or mental, arising out of persons incorrectly or inappropriately using or handling our products. We also assume no liability for any damage to an instrument or product from incorrect installation, handling or use of the product by the user.

Instructions

1. Using a sharp razor or scalpel, cut the 1/16" sample tube to desired length - The cut end must be square and burr-free. **Caution!** Be careful when handling and storing sharp objects.
2. Slide the cap and 1/16" ferrule over the cut end of the tube. Ensure that the 1/16" tube is protruding slightly above the ferrule face
3. Slide 1/8" cap nut over the nebulizer sample-inlet stem. Ensure that it is positioned as shown below.
4. Slip the 1/8" ferrule over the nebulizer sample-inlet stem. Note! Tapered portion of the ferrule must be pointing towards the 1/8" connector. Position the 1/8" ferrule with its face just below the protruding nebulizer stem.

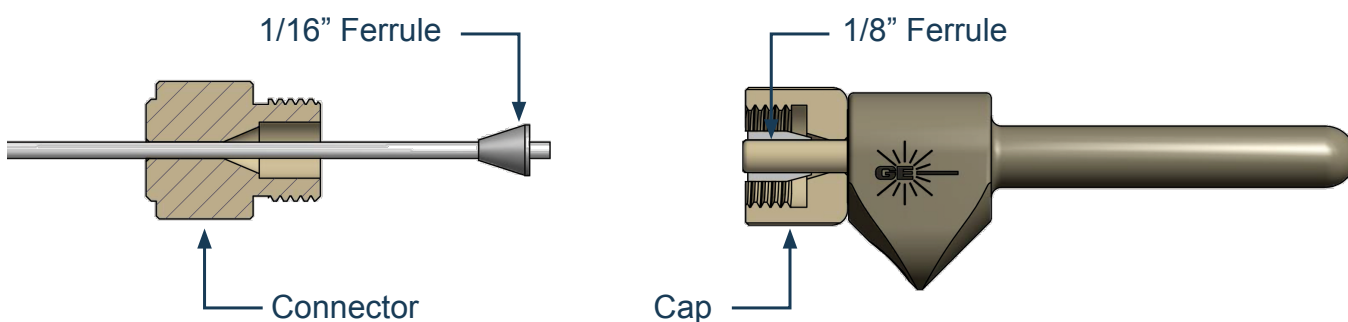

5. Bring together the cap (with 1/16" tube and ferrule) to the 1/8" connector (with Nebulizer and 1/8" ferrule). Hold cap and begin to screw together the connector until tight.
6. When finished, gently pull on the sample tube to ensure a firm connection has been made.

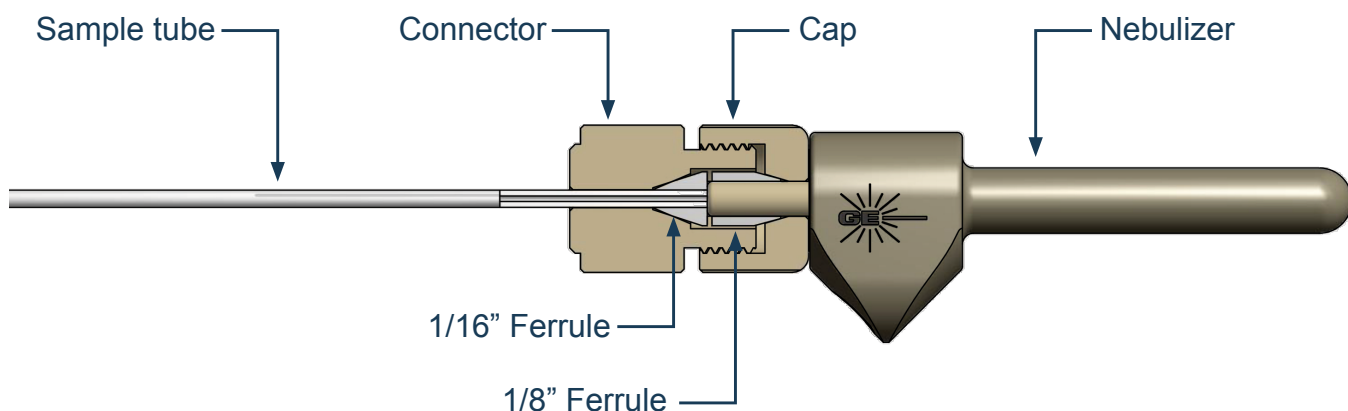

Caution! It is advised you wear safety gloves and glasses when handling glass.

Consumables

Replacement consumables can be purchased under the following part numbers.

Part Number	Description
70-803-1315	Tefzel Ferrule 1/16" (PKT 10)
70-803-1513	Tefzel Ferrule 1/8" (PKT 10)

For more information please visit www.geicp.com or contact one of our offices below.

Asia Pacific

6 Central Boulevard, Port Melbourne, VIC
3207, Australia

Telephone: +61 3 9320 1111
Email: enquiries@geicp.com

Americas

31 Jonathan Bourne Drive, Unit 7 Pocasset,
MA 02559, USA

Telephone: 508 563 1800
Email: geusa@geicp.com

Europe

Friedenbachstrasse 9, 35781 Weilburg,
Germany

Telephone: +49 6471 3778517
Email: gegmbh@geicp.com